

Standard Recommended 1x Dilution Scheme



Step 1

Take 25 ml bottle of Happy Cell[®] ASM 4X containing:


- 1X Culture Medium
- 4X Suspension Agent
- 1X Penicillin/Streptomycin



Step 2

Prepare 75ml bottle of Culture Medium containing:

- 1.35X FBS
- 1.35X Glutamine*
- 1X Penicillin/Streptomycin**

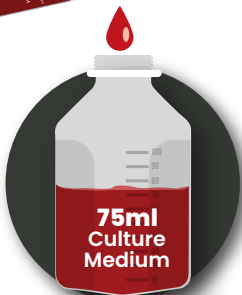

Step 3

Add the 25ml bottle of Happy Cell[®] ASM 4X to the 75ml bottle of culture medium. Gently mix the two together.

* At desired concentration ** If applicable